

# MAXPAR<sup>®</sup> Antibody Labeling Kit

## Lanthanide Labeling of Antibodies: *Pre-Load Method*



**CHEMICAL HAZARD:** Before handling any chemicals, refer to the Material Safety Data Sheet (MSDS) provided by the manufacturer, and observe all relevant precautions.

### Note:

- *This protocol has been optimized for a multitude of IgG isotypes and also works well for affinity purified polyclonal preparations. This protocol will NOT work with IgM antibodies.*
- *Each reaction is optimized for labeling 100 µg of antibody with a lanthanide metal.*

## Reagents and Materials

### Kit Contents (4 rxn):

- R-Buffer (6 mL)
- C-Buffer (5.5 mL)
- L-Buffer (1.4 mL)
- W-Buffer (8 mL)
- MAXPAR<sup>®</sup> Polymer (4 tubes)
- Lanthanide solution (20 µL)

### Kit Contents (40 rxn):

- R-Buffer (60 mL)
- C-Buffer (55 mL)
- L-Buffer (14 mL)
- W-Buffer (80 mL)
- MAXPAR<sup>®</sup> Polymer (40 tubes)
- Lanthanide solution (200 µL)

### Additional materials required:

- Antibodies to be labeled:  
Purified IgG or polyclonal: carrier-free (no BSA, Hydrolyzed protein, gelatin etc. for stabilization)
- Centrifugal Filter Unit: 3 kDa Amicon Ultra- 500 µL V bottom, Millipore Cat # UFC500396
- Centrifugal Filter Unit: 50 kDa Amicon Ultra- 500 µL V bottom, Millipore Cat # UFC505096
- Micro-centrifuge (*preferably 2 units to avoid overlap of certain steps requiring centrifugation*)
- Heat block Incubator or water bath (37°C)
- 0.5M TCEP: Bond-Breaker<sup>™</sup> TCEP solution, Pierce Cat#77720
- Aerosol Barrier (Filter) Pipette Tips
- PBS – based antibody stabilization solution: Antibody Stabilizer, Candor Biosciences Cat#131050, supplemented with 0.05% azide after purchase

### Storage:



- Buffers and Lanthanides: 4°C
- Polymers: -20°C with provided desiccant in a sealed container

### Important Notes Before Starting

- The MAXPAR<sup>®</sup> Polymer is moisture-sensitive. Equilibrate Polymer (stored at -20°C) to room temperature before opening to avoid moisture condensation.
- Use filter tips in all pipetting steps to prevent cross-contamination between metal stocks and reagents.
- Prior to starting the MAXPAR protocol, purified, carrier free antibody concentration should be verified by Nanodrop after blanking against **the buffer they are suspended in**. The composition of the buffer can be found on the technical data sheet supplied by the antibody vendor.
- Loading of the polymer and partial reduction of the antibody should be performed simultaneously (See Figure 1 at end of protocol). It is imperative, however, to not exceed recommended reduction time, or allow the partially reduced antibody to remain free of the loaded polymer.

Timing hh:mm (Approximate)	Outline of Protocol Steps
0:00	<b><i>Pre-Load the Polymer with Lanthanide</i></b>
	1. Spin polymer tube for 10 seconds in a microfuge to ensure the reagent is at the bottom of the tube
	2. Resuspend polymer with 95 $\mu$ L of L-Buffer
	3. Mix thoroughly by pipetting
	4. Add 5 $\mu$ L of Lanthanide metal solution to tube (final concentration: 2.5mM)
	5. Mix thoroughly by pipetting
	6. Incubate at 37°C for 30-40 minutes. During incubation, begin Buffer Exchange and Partial Reduction of Antibody
0:30	<b><i>Buffer Exchange and Partially Reduce the Antibody</i></b>
	7. Add 300 $\mu$ L R-Buffer to a 50kDa filter
	8. Add up to 100 $\mu$ g in up to 200 $\mu$ L of stock antibody to the R-Buffer in the filter (Note: if the stock antibody concentration is too dilute to add the desired amount of antibody, pre-concentrate it in the same filter prior to addition of the R-Buffer)
	9. Centrifuge 12,000 xg for 10 minutes at RT
	10. During centrifugation, dilute 0.5M TCEP stock to 4mM in R-Buffer by mixing 8 $\mu$ L of 0.5M TCEP stock with 992 $\mu$ L of R-Buffer. For each antibody being labeled 100 $\mu$ L of 4mM TCEP-R-Buffer is required.
	11. Discard column flow-through from centrifugation
	12. Add 100 $\mu$ L of the 4mM TCEP-R-Buffer to each antibody and mix by pipetting
	13. Incubate at 37°C for 30 minutes (do not exceed)
0:45	<b><i>Purify Lanthanide-loaded Polymer</i></b>
	14. Add 200 $\mu$ L of L-Buffer to a 3kDa filter
	15. Add the metal-loaded polymer mixture to the filter containing the L-Buffer
	16. Centrifuge 12,000 xg for 25 minutes at RT
	17. Add 300 $\mu$ L of C-Buffer to the filter and centrifuge 12,000 xg for 30 minutes at RT
1:15	<b><i>Purify the Partially Reduced Antibody</i></b>
	18. Retrieve the partially reduced antibody from the 37°C
	19. Add 300 $\mu$ L of C-Buffer to each 50 kDa filter
	20. Centrifuge 12,000 xg for 10 minutes at RT. A second microcentrifuge could be used at this step to avoid timing conflict with the polymer wash.
	21. Discard flow through
	22. Add 400 $\mu$ L of C-Buffer to the filter
	23. Centrifuge 12,000 xg for 10 minutes at RT
	<b><i>Retrieve the Partially Reduced Antibody and Lanthanide-loaded Polymer</i></b>
1:40	24. Retrieve 3 kDa filter containing the Lanthanide-loaded Polymer from the centrifuge and discard column flow through
1:55	25. Retrieve 50 kDa filter containing the Partially Reduced Antibody from the centrifuge and discard column flow through

Timing hh:mm (Approximate)	Outline of Protocol Steps
1:55	<b>Conjugate Antibody with Lanthanide-loaded Polymer</b>
	26. Using a pipette, re-suspend the lanthanide-loaded polymer in 60 $\mu$ L of C-Buffer
	27. Transfer the re-suspended contents to the corresponding partially reduced antibody in the 50 kDa filter
	28. Mix briefly by pipetting
	29. Incubate at 37°C for at least 60 minutes (up to 2 hours)
2:55	<b>Wash metal conjugated antibody</b>
	30. Add 300 $\mu$ L of W-Buffer to the antibody conjugation mixture
	31. Centrifuge 12,000 xg for 10 minutes
	32. Discard flow through
	33. Repeat 3 more times with 400 $\mu$ L of W-Buffer (for a total of 4 washes with W-Buffer)
3:35	<b>Recover Metal Conjugated Antibody</b>
	34. Add 50 $\mu$ L of W-buffer to the 50 kDa filter, pipette to mix and rinse the walls of the filter
	35. Invert the 50 kDa filter over to a new collection tube (see Figure 2)
	36. Centrifuge the inverted filter/collection tube assembly at 1000 xg for 2 minutes
	37. Remove the inverted filter from the collection tube, rinse walls of the filter with an additional 50 $\mu$ L of W-Buffer and replace it, inverted, back to the collection tube
	38. Centrifuge the inverted filter/collection tube assembly at 1000 xg for 2 minutes
3:40	<b>Yield Determination and Storage of Metal Conjugated Antibody</b>
	39. Quantify the conjugated antibody by measuring the absorbance at 280 nm against a W-Buffer blank (expected recovery is 60% )
	40. For storage, dilute the antibody to a final concentration of 0.5 mg/mL in a commercially available antibody stabilization buffer (supplemented with 0.05% sodium azide after purchase) and store at 4°C until ready to titrate
	41. Once the conjugated antibody has been titrated on the CyTOF it can be diluted to the optimum working concentration in stabilization buffer and stored at 4°C.

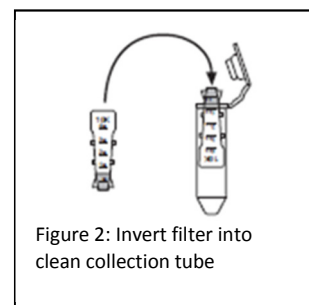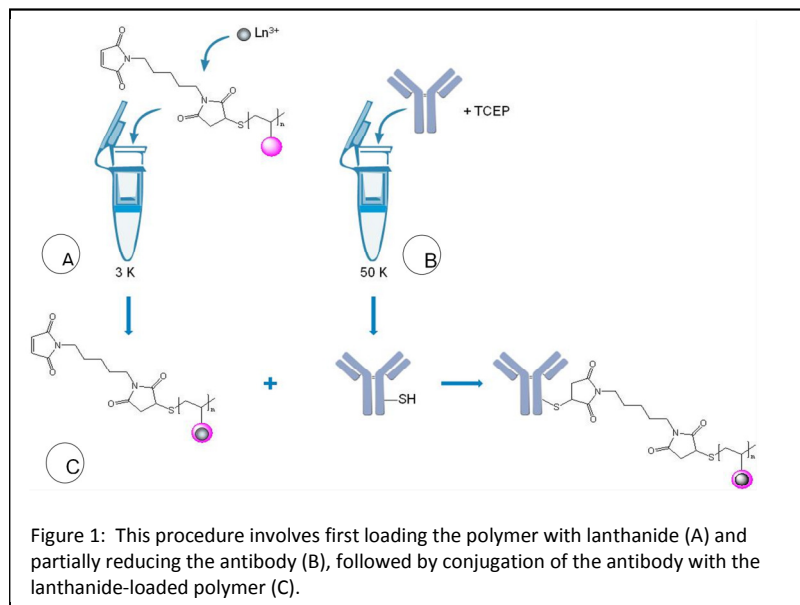

© 2012 DVS Sciences Inc. The trademarks mentioned herein are the property of DVS Sciences Inc. or their respective owners. For Research Use Only. Not for use in diagnostic procedures.

Information in this document is subject to change without notice. DVS Sciences Inc. assumes no responsibility for any errors that may appear in this document. This document is believed to be complete and accurate at the time of publication. In no event shall DVS Sciences Inc. be liable for incidental, special, multiple or consequential damage in connection with or arising from the use of this document.

DVS Sciences Inc. products are warranted to meet stated product specifications and to conform to label descriptions when used and stored properly. Unless otherwise stated, this warranty is limited to 6 months from date of sales for products used, handled and stored according to DVS Sciences' instructions. DVS Sciences' sole liability for the product is limited to replacement of the product or refund of the purchase price. DVS Sciences products may not be resold, modified for resale or used to manufacture commercial products without prior written approval from DVS Sciences Inc. DVS Sciences strives for complete customer satisfaction. If you are not satisfied with the performance of a DVS Sciences product, please contact DVS Sciences